

EloCheck 2.0 demo version

Please check this information before installing this software on your PC

- This free downloadable program package contains a demo version of the actual EloCheck 2.0 controller software from biotronix. The demo simulates in the background a virtual connected EloCheck unit so that almost all functions of the full software version are available for testing.
- System requirements: Microsoft NT, 98, 2000, ME or XP. For best performance a desktop resolution of 1024 x 768 Pixels (or higher) is required.
- Installation: Copy the self-extracting program 'EloCheck_demo.exe' to your computer and start it. The EloCheck software will be extracted and installed automatically. Click the 'EloCheck' button placed on your desktop after successful installation.
- To activate the connection to a virtual EloCheck unit press the 'Connect' button after starting the program.
- After pressing the 'Start' button (▶) the recording of a simulated growing culture starts. The whole simulation ends after about one hour. Restart the software to repeat the demonstration.
- Some examples of previously recorded data are free to download. For example you can find a recording of a 50 ml culture of *C. glutamicum* grown over six days within the file set 'Glutamicum1.exe'.
- For storing reports as *.pdf files the user need a pdf generating program on his computer. If no program as Adobe Acrobat, Adobe PDF etc. is present, the user can e.g. install 'FreePDFXP' / Ghostscript. Generation of pdf files can be done afterwards by selecting PDF-Maker as a virtual printer. By pressing 'print' PDF-Maker will open automatically.
- Please check the manual or the Help-menu for further details. If you need more information of if you have additional questions, do not hesitate to contact biotronix (info@biotronix.de)



↓ Connect

